

TEMPERATURE -20°C ~ +150°C





73 XI'AN LIB ENVIRONMENTAL SIMULATION INDUSTRY

Climatic Test Chamber Manufacturer And Supplier

Benchtop Temperature & Humidity Test Chamber

20%~98%RH



[Technical Parameter]

	Model	TH-50		TH-80
l	nternal Dimension (mm)	320*350*450		400*400*500
(Overall Dimension (mm)	800*1050*950		900*1100*1000
Interior Volume		50L		80L
Heat load		1000W		
	Temperature Range	A : -20 °C ~ +150 °C		
		B∶-40℃ ~ +150 ℃ C∶-70℃ ~ +150 ℃		
	Temperature Fluctuation	C: -/0℃ ~ +150 ℃ ± 0.5 ℃		
Parameter	Temperature Deviation	± 2.0 ℃		
mete	Humidity Range	20% ~ 98% RH		
Ч.	Humidity Deviation	± 2.5% RH		
	Cooling Rate		1 ℃	/ min
	Heating Rate		3 °C	/ min
		Cooling system Mechanical compression refrigeration system		
	Cooling	Refrigerating unit French TECUMSEH compressor		
		Refrigerant		
	Heating Element	Nichrome heater		
		Programmable color LCD touch screen controller		
	Controller	Ethernet connection, PC Link		
Stru		Humidifier	External isolation, s	tainless steel surface evaporation humidifier
tructure	Humidity	Dehumidification		Evaporator
Ire	Concer	Temperature Sensor	PTR Platinum Resist	tance PT100Ω/MV A-class, accuracy 0.001 $^\circ\!\!\!\!\!\!^\circ$
	Sensor	Humidity Sensor		Dry and wet bulb sensor
	Build-in Water Tank(mm)	150*350*180		
	View Window Size(mm)	120*150		
	Air Circulation	Centrifugal wind fan		
	Safety Device	Humidifier Dry-combustion protection; Over-temperature Protection; Over-current Protection;		
	Salety Device	Refrigerant High-pressure protection; Water Shortage Protection; Earth leakage Protection		
	Exterior Material	Steel Plate with protective coating		
Material	Interior Material	SUS304 stainless steel		
erial	Thermal Insulation	Polyurethane foam and insulation cotton		
	Observation Window	Interior lighting, double-layer thermo stability silicone rubber sealing		
	Standard Configuration	1 Cable hole (Φ 50,) with plug; 2 shelves		
	Power Supply	220V 50Hz / 380V 50Hz		
Maximum Noise		65 dBA		
Environmental Conditional		5℃~+35 ℃ ≤85% RH		
 Low temperature can con 		l down to -86℃. □ Standard & customized		

 $\hfill\square$ Low temperature can cool down to -86 $^\circ\!\mathbb{C}.$

Standard & customized

- The Refrigerant can be replaced by R449A/R508B.
- □ The heat load can absorb 1000W in -20°C.

- * pay attention to the environmental condition during installing






1.Workroom

The internal material is 304 stainless steel, mirror surface, rust-proof to high and low temperature and moisture corrosion;

Nice ruggedness and long life.

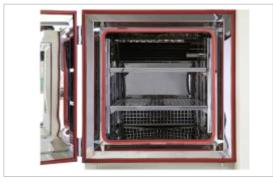

2.Sample Shelf

SUS304 stainless steel punching sample shelf 2pcs; Standard load of 50KGS for each layer of sample shelf; The height of the shelf is adjustable and can be removed directly to facilitate the various sizes of test samples. Conducive to the air circulation.



3.Cable Hole

A standard test hole, located at the left side of the work room, is provided with a sleeve inside to prevent moisture from entering the insulation layer. Equipped with flange. The standard inner hole diameter is 50mm.

4.Viewing Window

Double layer insulating glass 8cm thick, made of tempered glass.

The conductive film is located on the interior glass to prevent window frosting, built-in LED light for the work room lighting ;can clearly observe samples.

The light control button is located below the controller.



5.Controller

PID programmable color touch screen controller, network connection computer.

Can create 120 programs, 100 segments.

LIB also can preset program into the controller based on user testing requirements.

The set system language is English for standard.

HRIN PROG PTOO1 HEXT	HAIN PROG PTOO1 NEXT
TEMP * -20.00 °C	TEMP * 85.00 °C
HUMI - %	BP 85.0 85.0 %
PROG/SEG: 001/002 RUN TIME: 00H02M 19-18-09 10 59 RH JUMP HOLD STOP	PROG/SEG: 001/002 RLN TIME: 00H02 19-10-09 10 59 RH JUHP HOLD STOP

6.Build-in Water Tank

The water stored in the built-in water tank meets the needs of the test. When there is a low water level, it will alarm to indicate that water needs to be added. At this time, it is necessary to add enough water till to the high water level.

7.Castor

Install 4 castors for ease moving, and with brakes function. Caster height adjustable.



8.Insulation

10cm thick polyurethane foam and insulation cotton; Better insulation performance, degradation resistance, environmental protection, noise reduction.



9.External Materials

A3 steel plate with galvanized coating; Electrostatic treatment; High and low temperature corrosion resistance; High hardness, anti - impact Very high safety factor; Color can be customized.


Soft silicone plug seal; Metal enclosure closed, beautiful designed.





Structures

Core Function

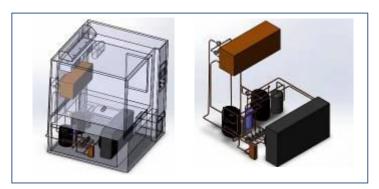

Cooling System


After the liquid refrigerant absorbs the heat of the cooled object in the evaporator, it is vaporized into low-temperature and low-pressure steam, sucked by the compressor, compressed into high-pressure high-temperature steam, discharged into the condenser, and radiated to the cooling medium in the condenser, condensed into a high-pressure liquid, throttling through a throttle valve into a low-pressure low-temperature refrigerant, and again entering the evaporator to absorb heat and vaporize, to achieve the purpose of cycle refrigeration.

In this way, the refrigerant completes a refrigeration cycle through four basic processes of evaporation, compression, condensation, and throttling in the system.




Humidity System

When the current humidity is less than the set point humidity, the equipment is automatically humidified. The humidification of is automatically heated by the humidification tank. When the water is heated, steam is generated, and then the steam is injected into the working room to increase the test humidity.

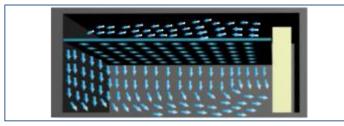

Air Circulation

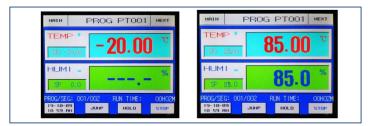

Controller System

uniformly distributed through the air outlet. Air circulation adopts air outlet at top and air return at bottom, and the wind speed and pressure are in compliance with the test standard, and the temperature is stabilized at the moment of switching. The high-quality centrifugal fan is used to strongly supply air circulation, making the temperature distribution in the test area uniform.

The centrifugal fan is installed at the rear of the chamber body, and the air is

The PID controller as the main control unit to command, operate, detect and redistribute the various components of the equipment to achieve maximum effective use.

The temperature control adopts P.I.D. S.S.R. system synchronous coordinated control, which can improve the stability and life of the control components and interface.

Screen display function: LCD display, which can display test conditions (including temperature section, cycle number, running time and remaining time, etc.).



Temperature and Humidity Sensor

PT-100 Class A sensor, real-time accurate detection and display of temperature changes at 0.001 degrees.

A wet gauze and a real-time temperature, converted by temperature difference, showing real-time relative humidity.

Ver Temperature Drotestian

Over Current Protection

lumidifier Dr. Cemphustien Dretestien

Water Shortage Protection

Earth Leakage Protection Phase Sequence Protection Refrigerant High Pressure Protection

Protection System

Sample Shelf	 Quantity can be increased Special sample shelf can be customized
Cable Wire	 2.5m 3m
Test Hole	 Diameter 100mm Diameter 125mm Shape and size can be customized
Capacitive Temperature And Humidity Sensor	VAISALAROTRONIC
Observation Window	 Size can be customized Inner glass door Hand-in ports
Reverse Osmosis Unit	
Controller	 WiFi connection Mobile APP control Multi-language controller interface: Russian, French, Polish, Romanian, German

Shipping & Packaging

Environment Conditions


Temperature: +5°C ~ +35°C Relative humidity: ≤85%; Pressure: 86 KPa—106 KPa

Space Conditions

Sufficient space for ventilation



Install Condition

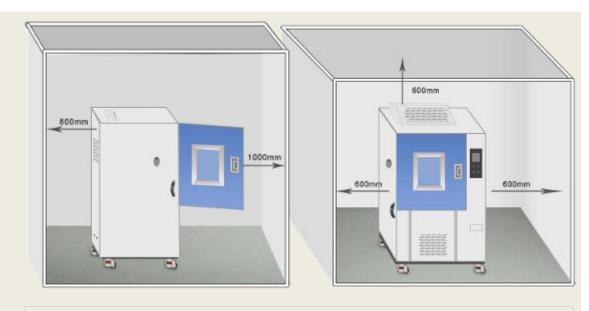

*Safety Instruction: 1.Prohibited to test explosive, inflammable and high corrosive substance 2.Chemical exposure to the equipment is prohibited 3.Equipment must be safety on the ground to avoid electrostatic induction

Packaging



First, seal chamber with waterproof plastic film. Protect chamber from seawater corrosion. Second, buffer foam is placed in the four corners of the chamber. It is used for fixing equipment to prevent shaking and damaging chamber during transportation.



Plywood: standard wood export packaging. The wooden box is fixed by sheet metal to prevent damage during transportation.

Gross weight: 180kgs~ Package size: 1170*960*1150 mm

Warranty Condition

Our company will repair the product, if the product, the material of the parts, the design and manufacturing of the products raised hardware problems caused by product itself rather than human error within three years warranty period since the date of dispatching by the customer. Our company repairs the products, but will collect the basic costs of the spare parts after the warranty period, but service is free always.

Shipping



Material: Export standard wooden box *Can be used for Sea, Air, Railway, Truck and multimodal transport.

Warranty & Service

- 3 Years Warranty, Lifelong Follow-up Services

Professional After-sales Team If you have any questions during operation, we will supply solution to you with 24 hours.

How to Service

1. At first, our test chambers are produced based on 20 years product lifetime.

Normally once test chambers have problems, we judge the problems, and send spare parts to our customers, and teach them how to change new parts on by email or video, all spare parts and shipping cost (by DHL, TNT, and FedEx) paid by LIB.

2. f the customer needs our engineer on-site service, they only need pay the ticket accommodation to our Engineers, service is for free.

3. If products still can't use after our engineers repair, we will produce a new test chamber (same as the old one) to our customers with no charge.



PROVIDING TEST SOLUTION, PROVIDING TEST EQUIPMENT

Xi'an LIB Environmental Simulation Industry is a lead provider of environmental test chamber in China, with its own brand (LIB) design, production, sales and service since 2009. We continually update technology and develop new products for customer's needs.

Our main products included temperature and climate test chambers, corrosion chambers, weathering testers, IP dust and rain chambers, ozone test chamber, noxious gas SO2 H2S chamber, walk-in chambers. We provide test chamber, we provide test solutions. Standard and customized products to meet different customer needs.


By 2018, global market has spread to 51 countries to USA, Canada, Mexico, Brazil, Peru, Russia, German,UK,France,Finland,Netherlands, Poland,Switzerland,Thailand,Philippines, Singapore, Malaysia, Australian, South Africa etc. and the market continues to expand. At present 7 Tier-one agents around the world provide LIB products, installation and maintenance service for customers. Make things simple and convenient.

Contact Us

PARTNER

Website: www.lib-industry.com Add: No.5 Zhangba First St, Xian High-Tech Area, Shaanxi ProvinceP.R.China 710065 Tel: +86 29 68918976 WhatsApp: +86 18700875368 Skype: serena_lib Email: serena@ lib-industry.

Sales Team: <u>sales@lib-industry.com</u> After Sales Team: asd@lib-industry.com Shipping Team: vip@lib-industry.com Finance Team: account@lib-industry.com

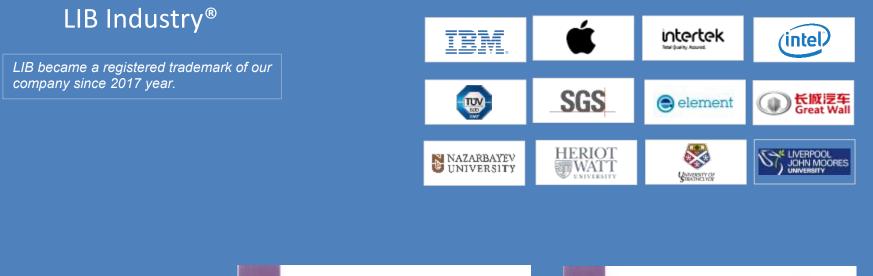







Contificate of Conformity

ROHS

Ξ	Applicant	: 3
C2	Menulacturer	: 3
	Trademark	1.6
	Product	1.1
	REN	: 7
Ĩ		1 1 1
0		Ť
0	Test Standard	
NFORMI	The cartificate mentioned pro possible to use 20084/2EC as report member:	ducts • CE d LVI
TY	6/F, Build Shenz E-mail	thei.

		idian Industry High-Tech Area, Shaamii Province,
		ntion Industry High-Tech Anes, Shaamsi Province,
1.18		
Temperatur	e Humidity Test Cha	mber
TH-1000, T TH-5000, T TH-10000, T T-1004, T-1	H-100C, TH-225A, TH H-500C, TH-600A, TH TH-1000C, T-60A, T-5 BOIL, T-100C, T-225A,	A. TH-865, TH-86C, TH-160A, 4-2558, TH-255C, TH-608A, 4-8508, TH-800C, TH-608A, 4-8568, T-80C, T-408A, T-60C, 17-2258, T-30A, T-608A, 18, T-826C, T-608A, 18, T-826C, T-1030A, T-1030B,
EN ISD 121 EN 60204-1	00:2010 :2006+A1:2009+AC:5	2010
ts. If doars in 5 marking to	of imply an assessme demonstrate the co 2014/35/EU. It is onl	union of a sample of the above ent of the whole production. If is impliance with this MD Directive by valid is connection with the test
E		(ALA) Manager Jon 272 (\$13
3, Zhoulong 1 , Guangdon;	, China. Tel:4000-83	Go., Ltd. an Street, Longgang Diatrict, 76-382 0766-32925541 h; http://www.bko-kab.com

Applicant	No.5 Zhang	Siwironmontal Birmlat gba First Street, Xi'an H PR China 710085	ion Industry gh-Tech Acea, Shaaree
Manufacturer	No.5 Zhang	Invironmental Simulat pa First Street, Xi'an Hi 'R China 710055	
Trademark	3.10		
Product	: Temperatu	ne Humidity Test Char	sher
56.74	TH-1008, 1 TH-5008, 1 TH-10008, T-80C, T-9	CH-109C, TH-225A, TH- CH-509C, TH-800A, TH- TH-1000C, T-60A, T-50 80A, T-100E, T-500A, T-50 80A, T-100E, T-500C, T-500A	, TH-808, TH-80C, TH-100 2258, TH-225C, TH-560A, 6008, TH-800C, TH-560A, 8, T-50C, T-80A, T-808, -255A, T-2258, T-235C, , T-8008, T-800C, T-1080A,
	a second a		
The EUT desc		been consolidated by u	
he EUT dasa he council I o demonstrati	ribed above has toHS directive the compliance	been consolidated by u 2011/65/EU. If is p	ossible to use CE mark
The EUT dasa the council I to demonstrate	ribod above has toHS directive r the compliance port number: BK	been consolidated by u 2011/85/EU. If is p with this RoHS Deech	s and found is compliance to contribut to use CE strack, w. E is only wake in connect
The EUT dosc the council J to demonstrate with the test is	ribod above has toHS directive r the compliance port number: BK	been consolidated by u 2011/85/EU. If is p with this RoHS Deech	osaible to use CE mark